

ExCell Bio

ResiQuant® CHO HCP ELISA Kit-FAST

For Research and Manufacturing Use

Not Intended for Diagnostic and Therapeutic Use

User Manual

Catalog Number CRH00-3051S

CRH00-3051

CRH00-3052



ExCell

PRODUCT DESCRIPTION

The expression of therapeutic proteins in the CHO cell lines is economical and convenient for commercial drug

production. However, the host cell proteins (HCPs) from CHO cells as a type of concomitant impurities are one

of the major concerns in terms of drug safety. These impurities may diminish the efficacy of drug therapy and

cause side effects or unexpected immune reactions. Therefore, control and monitor of the residual CHO HCP

impurities is world widely required by all corresponding regulatory authorities. This product is a sandwich

ELISA kit. The antibodies were generated from rabbits and goats against the antigens harvested from the

fermentation supernatant of CHO cells, cultured in the medium free of proteins. This kit can be used throughout

the manufacturing process and for product release.

For the first time, it is recommended to conduct a suitability study to confirm the sample matrix effect and

determine the appropriate dilution conditions for testing.

PRINCIPLES OF THE TEST

The assay employs a sandwich immunoassay methodology for the quantitation of CHO HCP. Microplates are

coated with antibodies that exhibit specificity for CHO HCP. Anti-CHO HCP antibodies labeled by HRP and

standards or samples are added to the wells simultaneously. During incubation, CHO HCP in standard and

samples forms immune complexes with the antibodies. After washing away unbound substances, the

chromogenic substrate 3,3',5,5'-tetramethylbenzidine (TMB) is introduced, initiating an enzymatic reaction. This

reaction involves HRP-mediated oxidation of TMB in the presence of hydrogen peroxide (H2O2), yielding a

blue-colored product that exhibits maximum absorbance at 655 nm. The enzymatic reaction is subsequently

halted by the addition of a stop solution, which turns the color to yellow with peak absorbance at 450 nm. The

optical density (OD) value at 450 nm, measured by an enzyme-linked immunosorbent assay (ELISA) reader, is

proportional to the concentration of CHO HCP within the samples. Quantitative determination of CHO HCP is


achieved with a simultaneously prepared standard curve, using the standards supplied within the kit.

Web: www.excellbio.com Tel: +86-021-3367 6611 Email: marketing@excellbio.com

1



Schematic diagram:

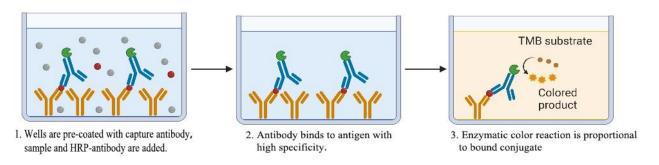

PRODUCT PERFORMANCE

- 1. Sensitivity: The limit of detection (LOD) for CHO HCP is 0.5 ng/mL, and the low limit of quantification (LLOQ) is 1.0 ng/mL.
- 2. Precision: The coefficient of variation (CV) for sample concentrations within the quantitation range is no more than 20% for both intra- and inter-assay.
- 3. Specificity: No cross-reactivity has been observed with E. coli, Vero, HEK293 whole cell lysates.

| PRODUCT APPLICATION

This product is a general assay kit designed for the quantitative detection of CHO HCP impurities in samples.

PRODUCT SPECIFICATIONS

Catalog Number	Name	Size
CRH00-3051S	ResiQuant® CHO HCP ELISA Kit-FAST	48T
CRH00-3051	ResiQuant® CHO HCP ELISA Kit-FAST	48T
CRH00-3052	ResiQuant® CHO HCP ELISA Kit-FAST	96T



REAGENTS PROVIDED

Name	96 Tests	48 Tests	Store				
CHO HCP Standard							
Standards at 250, 100, 40, 16, 6.4, 2.6, 1	8	8	2°C to 8°C				
and 0 ng/mL, 1 mL/vial							
CHO HCP 2G Microplate	8× 12	8× 6	2°C to 8°C				
100× CHO HCP HRP-Antibody	60 μL	30 μL	2°C to 8°C				
CHO HCP Assay Diluent	2× 25 mL	25 mL	2°C to 8°C				
20× Wash Buffer Concentrate	30 mL	30 mL	2°C to 8°C				
Substrate Solution	12 mL	6 mL	2°C to 8°C (Light-Sensitive)				
Stop Solution	12 mL	12 mL	2°C to 8°C				
Plate Sealer	3	2	/				

INSTRUCTION FOR USE

I. Materials required but not provided

- 1. Microplate reader capable of measuring absorbance at 450 nm (If wavelength correction is needed, additional wavelength at 630 nm or 570 nm is required).
- 2. $0.5-10 \mu L$, $2-20 \mu L$, $20-200 \mu L$, $100-1000 \mu L$ adjustable micropipettes with disposable tips.
- 3. Microplate shaker.
- 4. Deionized water.

II. Sample collection

- 1. Samples should be clarified and any precipitate should be removed by centrifugation.
- 2. The sample should be diluted appropriately with the CHO HCP assay diluent according to the pre-determined condition (It is recommended to complete the applicability study of the kit to confirm the appropriate dilution conditions for sample detection).

III. Preparation of reagents

- 1. It is recommended to remove the assay kit from the refrigerator 20 minutes in advance to allow it to equilibrate to room temperature.
- 2. Dilute 20× Wash Buffer Concentrate with deionized or distilled water to prepare the wash buffer. Store the

ExCell

unused solution at 2°C to 8°C.

3. CHO HCP HRP-Antibody Working Solution: Freshly prepare every time the working solution by diluting

100× CHO HCP HRP-Antibody with CHO HCP Assay Diluent to 1× in an amount sufficient advance 30

minutes for the current experiment.

IV. Washing

1. Hand-wash: add 300 µL of wash buffer into each well, allow it to sit for 10 seconds, then pour off the liquid

in the well. Invert the plate and blot it against clean paper towels. Wash the plate 5 times.

2. Automatic washing machine: set the volume to 300 µL, and the interval between aspiration and dispensing is

10 seconds. Wash the plate 5 times.

V. Assay procedure

1. Bring all reagents and samples to room temperature before use. Remove excess microplate strips from the

plate frame, return them to the foil pouch containing a desiccant pack, reseal tightly and store at 2°C to 8°C.

2. Set the blank well (If the plate is measured by dual-wavelength, the blank may not be necessary).

3. Prepare samples and standards in advance.

4. Add 100 μL different concentrations of standards or samples to the microplate wells respectively, then add

50 μL per well of CHO HCP HRP-Antibody Working Solution immediately. Seal the wells with the plate

sealer. Incubate at 37°C and shake for 120 minutes, using a microplate shaker (500 rpm).

5. Remove the liquid from the wells and wash the plate 5 times.

6. Add 100 μL of Substrate Solution to the wells, Incubate at 25°C in the dark for 15 minutes.

7. Add 100 µL of Stop Solution to the wells. Read the optical density (OD) of each well within 10 minutes

using a microplate reader set to 450 nm. If wavelength correction is available, the OD value at additional

wavelength (630 nm or 570 nm) should be collected. For single wavelength mode, the OD value of blank

well shall be subtracted from that of each standard or sample.

VI. Assay Procedure Summary



Add 100 μ L standards or samples to each well, then add 50 μ L CHO HCP HRP-Antibody Working Solution immediately. Incubate 120 minutes at 37°C, 500 rpm.



Remove the liquid and wash the plate 5 times.



Add 100 μL Substrate Solution to each well. Incubate at 25°C in the dark for 15 minutes.



Add 100 µL Stop Solution to each well. Read at 450 nm within 10 minutes.

| ANALYSES OF RESULTS



- 1. Create a standard curve by plotting the mean optical density (OD) of each standard concentration as the y-axis against the standard concentration on the x-axis. It is recommended to use a four-parameter fitting equation for the standard curve. Calculate the concentration of samples based on their OD values.
- 2. The coefficient of determination (R²) for the standard curve should be no less than 0.99. Discard those abnormal data, the deviation between the back calculated and the theoretical concentration of each standard should be within ±20% (±25% for the upper or lower limits of quantitation).
- 3. If the OD value of the sample exceeds the high end of the standard curve, the sample should be appropriately diluted before further measurement. Be aware that the dilution factor should be counted with back calculated concentration together.
- 4. The concentration of samples should be calculated with the standard curve in the parallel assay.

Representative Standard Curve:

Standard	OD450-630			Concentration (ng/mL)			3						
(ng/mL)	1	2	Average	1	2	Average	2.5						
250	2.606	2.493	2.550	261.0	239.2	250.1]						
100	1.515	1.465	1.490	102.9	97.8	100.3	2						
40	0.780	0.770	0.775	40.0	39.3	39.6	05.						
16	0.407	0.399	0.403	16.3	15.9	16.1	1.5		/				
6.4	0.217	0.231	0.224	6.2	6.9	6.6	1						
2.6	0.140	0.143	0.142	2.6	2.7	2.6		1					
1.0	0.104	0.104	0.104	1.0	1.0	1.0	0.5						
0	0.076	0.074	0.075					*					
\mathbb{R}^2	0.99999						0	0	50 Co	100 oncentration	150 (ng/mL)	200	250

Note: This graph is provided for reference purpose only. Data analysis was using ELISA Calc software with a four-parameter fitting model.

NOTES

1. Store the reagents at 2° C to 8° C.

ExCell

2. The concentrated HRP-Antibody is supplied in a small volume. Turbulence or inversion happened during

transportation may cause spray of the solution elsewhere on the wall or the cap. Therefore, please spin the

tubes shortly before use to harvest all liquid to the bottom of the tube.

3. In case crystal sediment appears in the 20× Wash Buffer Concentrate, warm it up at 37°C to dissolve the

sediment before use.

4. Do not mix or substitute reagents with those from other lots or different kits.

5. Ensure thoroughly mixing in preparing the solutions to ensure the reactions are consistent.

6. It is recommended to test all samples and standards in duplicate form at least.

DISCLAIMER

1. The product should be used according to the instructions in the manual. If the experimenter fails to operate

according to the instructions, our company will not be responsible for any deviation in product performance

caused by this.

2. The product is only used for scientific research and commercial production, and is not suitable for clinical

diagnosis and treatment. Otherwise, all consequences arising shall be borne by the experimenter, and our

company shall not be responsible.

Web: www.excellbio.com Tel: +86-021-3367 6611 Email: marketing@excellbio.com

7